


Instructions for Accessing Network Version WinList 3D

(a 40 reusable user license network share for Flow Cytometric analysis)

Click on :START” lower left of your desktop


In the box displayed after selecting Run, type [\\flowcyto\flowresearch\\$\VeritySoftwareHouse](\\flowcyto\flowresearch$\VeritySoftwareHouse) and select “OK”.

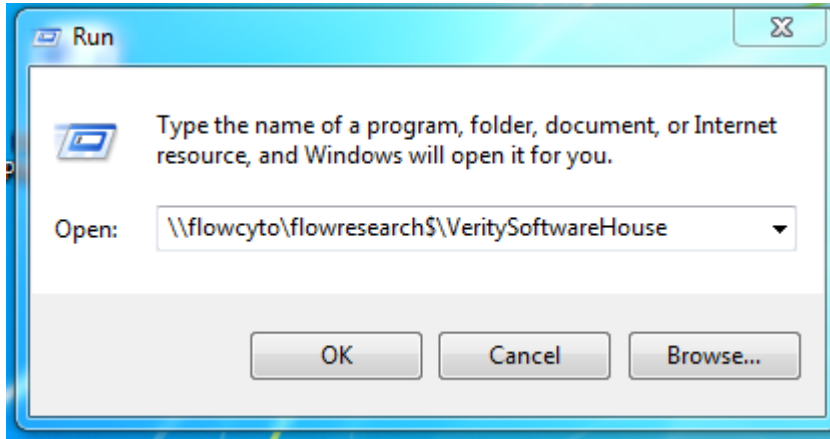

Note: You must be a member of this folder group to gain access.


Please E-mail ed.podniesinski@roswellpark.org to be added to the group.

This is the shared folder location where WinList3D_NetworkSetup resides.

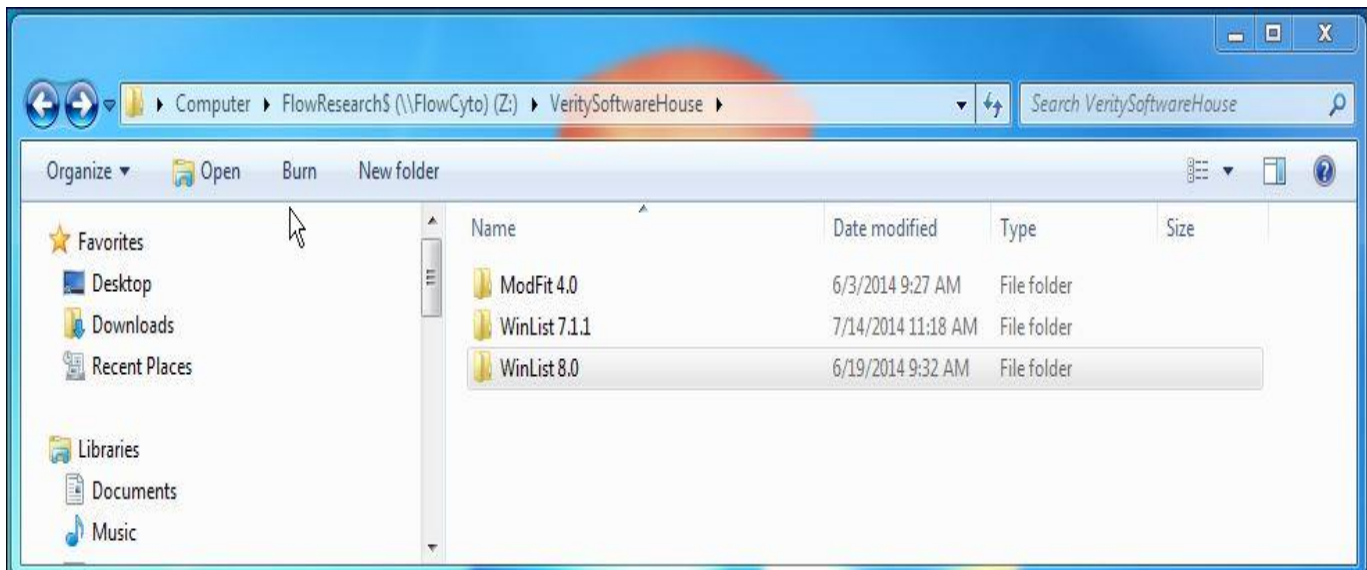

Open WinList 3D 7.1.1 or WinList 8.0 Network folder by double clicking on it:

Here's what you should see for WinList 7.1.1 directory contents:

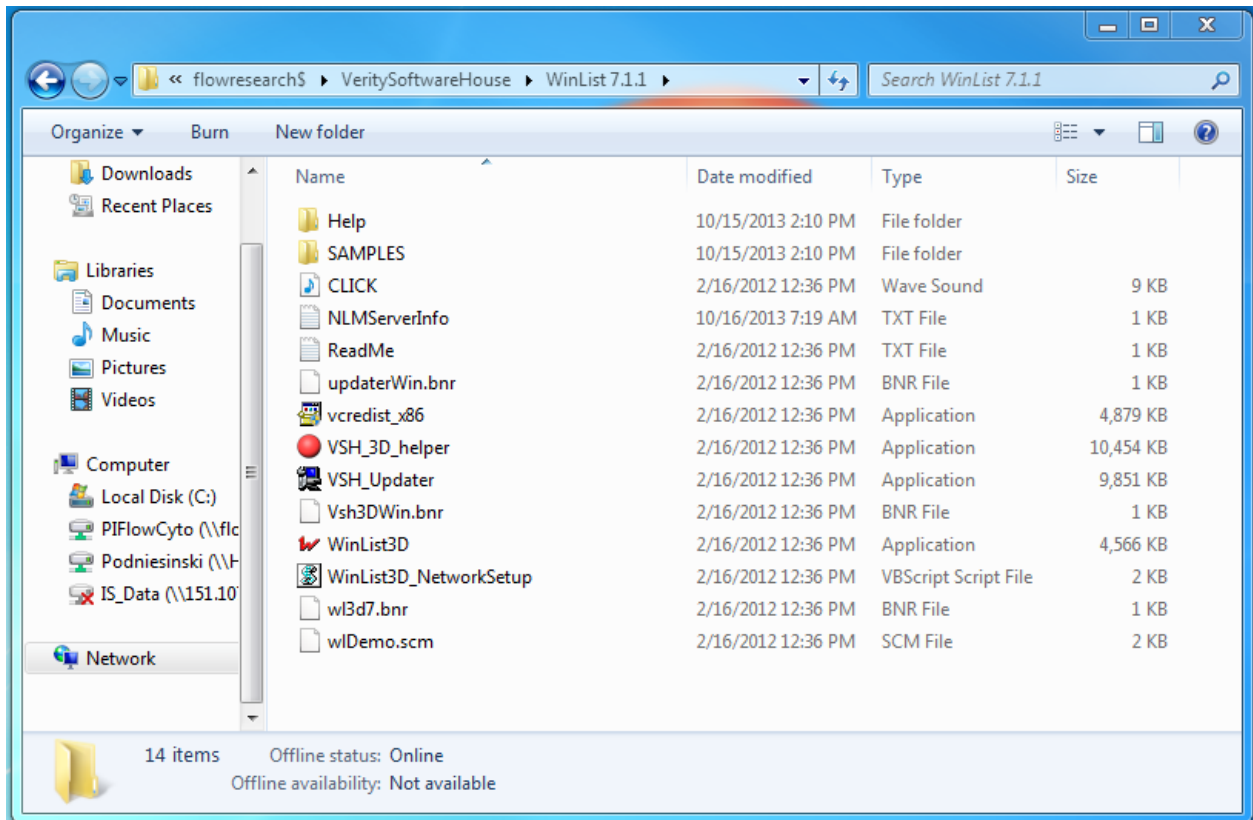

Right click on WinList3D.exe and within the popup menu, select "Send to" Desktop (Create Shortcut).

This will create a shortcut on your Desktop. You may want to rename it to WinList 7.1 Network.

For WinList 8.0 temporary Admin rights are needed to run the WinList3D_NetworkSetup script file. Once run once, computer Admin rights are not needed.

Running this script will create a desktop WinList icon automatically.

WinList 8.0 directory view, launch the highlighted automated script by double clicking:

WinList3D	3/18/2014 11:07 AM	Application	7,753 KB
WinList3D_NetworkSetup	3/17/2014 11:19 AM	VBScript Script File	2 KB
wl3d8.bnr	2/10/2014 3:56 PM	BNR File	1 KB

Notes:

WinList 8.0 version must use WinList3D_NetworkSetup script file but this requires temporary Admin rights and only needs to be run once.

WinList 7.1.1 Usually, using the first set of instructions in creating the Desktop shortcut works sufficiently without obtaining Temp Admin rights.